

Task

To examine the working principle of a potentiometer by means of a model of a potentiometer and subsequently to vary the brightness of a filament lamp by use of a commercially available potentiometer.

Equipment


| | | |
|---|----------|-----|
| Plug-in board | 06033.00 | 1 |
| Wire building block | 39120.00 | 4 |
| Lamp holder E10 | 17049.00 | 1 |
| Potentiometer, 250 Ω | 39103.21 | 1 |
| Universal holder | 39115.02 | 2 |
| Connecting cable, 50 cm, red | 07314.01 | 2 |
| Connecting cable, 50 cm, blue | 07314.04 | 2 |
| Crocodile clips, bare, 1 from 10 | 07274.03 | (1) |
| Filament lamp, 4V/0.04 A, E10, 1 pc. | 06154.03 | (1) |
| Constantan wire, d = 0.2 mm, need approx.. 30 cm | 06100.00 | (1) |
| Multi-range meter | 07028.01 | 1 |
| Power supply, 0...12 V-, 6 V~, 12 V~ | 13505.93 | 1 |
| Ruler | | |

Set-Up and Procedure

First experiment

- Connect up the circuit as shown in Fig. 1; fix the constantan wire so between the universal holders that it does not sag.
- Connect the crocodile clip on the freely movable wire connected to the voltmeter to the universal holder on the right.
- Select the 1 V- measurement range.
- Set the power supply to 0 V, then switch it on.
- Carefully increase the power supply voltage until the voltmeter shows 1 V.
- Measure the length l of the inserted piece of wire and note the measured value in Table 1.
- Move the crocodile clip to connect the voltmeter successively to different positions on the wire (e.g. about $\frac{3}{4}$, $\frac{1}{2}$, $\frac{1}{4}$ along the length of the wire), measure at each position the length of wire held and the voltage across this length of wire. Note the measured values for l and U in Table 1.
- Set the power supply back to 0 V and switch it off.

Fig. 1

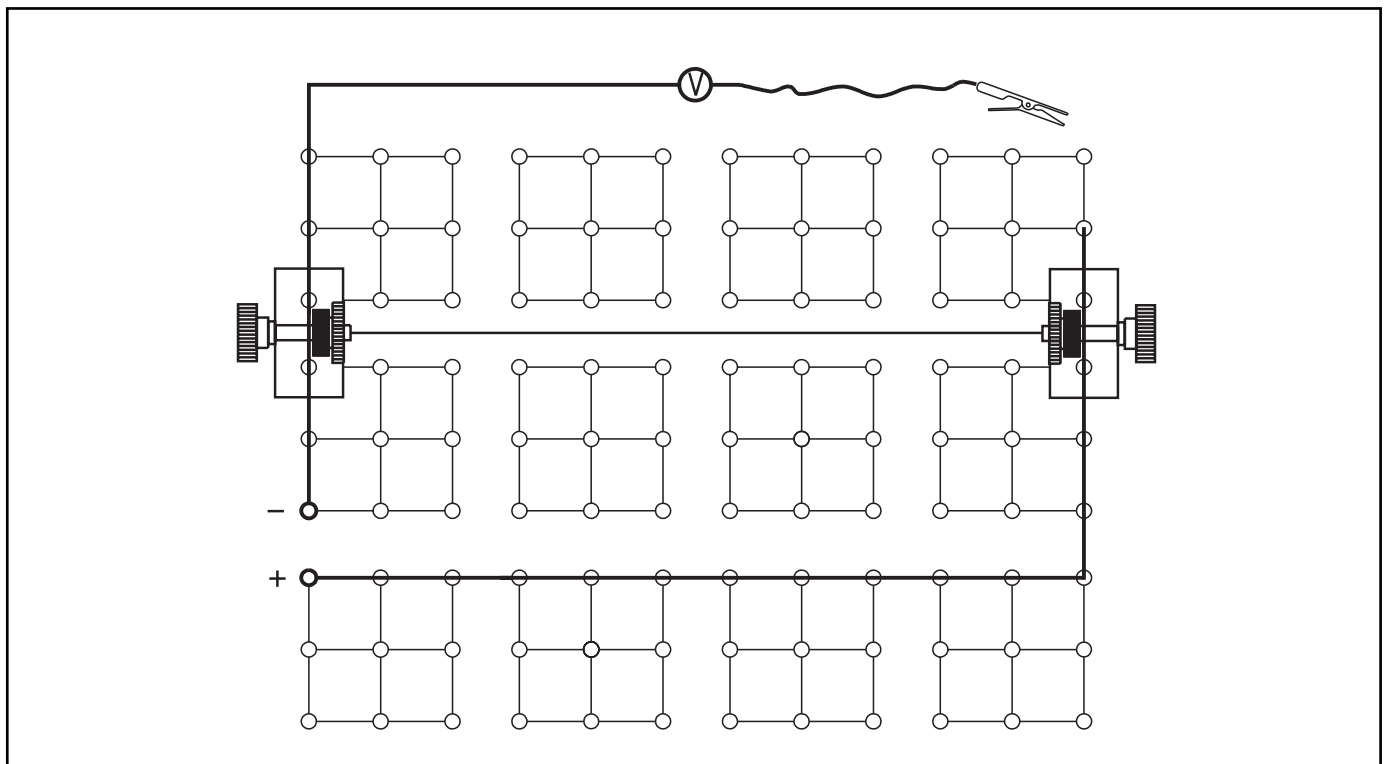

Second experiment

- Set up the experiment as shown in Fig.2; set the rotary knob of the potentiometer to 0; do not insert the filament lamp at first.
- Select the 10 V- measurement range.
- Switch on the power supply and set it to about 4 V.
- Slowly turn the potentiometer knob fully to the end stop and then back to 0; observe the deflection of the voltmeter while doing this.
- Note what you observed under (1).
- Complete the circuit by fitting in the filament lamp.
- Slowly turn the potentiometer knob from the 0 mark to the 10 mark and then back to 0, and observe the lamp while doing so.
- Note what you observed under (1).
- Set the power supply to 0, then switch it off.

(1)

[illegible]


(2)

Observations and Measurement Results

Table 1

| $\frac{I}{m}$ | $\frac{U}{V}$ | $\frac{U/I}{V/m}$ |
|---------------|---------------|-------------------|
| | 1.00 | |
| | | |
| | | |
| | | |

Fig. 2

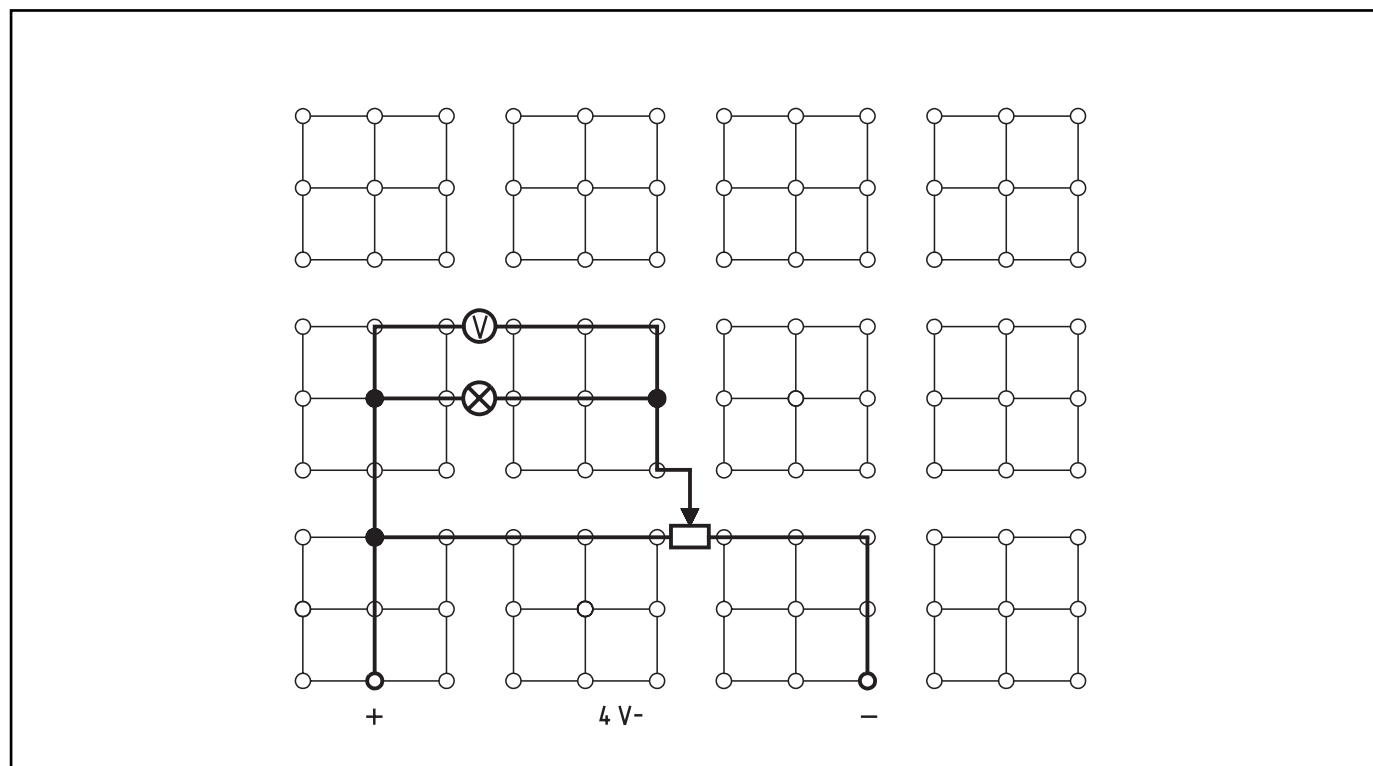



Evaluation

1. Plot a graph showing the dependence of the voltage U on the length of wire held l (Fig. 3). Which relationship between U and l can you derive from Fig. 3?

2. Form the quotient U/l for each pair of measured values obtained and enter these in column 3 of Table 1. Formulate the relationship between U and l mathematically.

Fig. 3

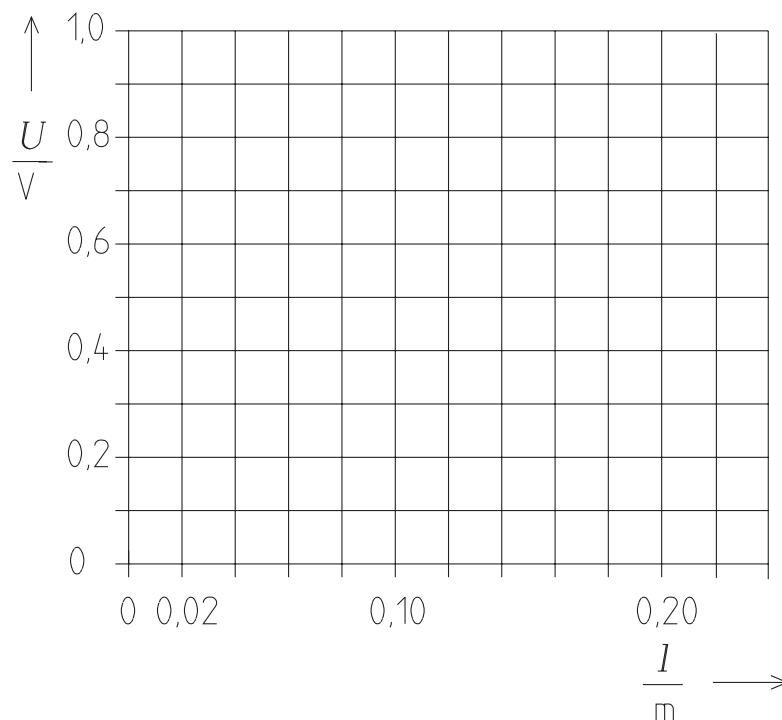

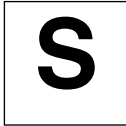

EEP
2.9

How does a potentiometer function?



3. The function of a potentiometer is given from the relationship determined under 2 and the observations noted under (1) and (2). Describe what a potentiometer can be used for.

.....

.....

.....

.....

.....

.....

.....

.....

.....

.....

.....

.....

.....

.....

.....

.....

.....

.....

.....

.....

.....

4. Give examples of where a potentiometer is used in practice.

.....

.....

.....

.....

.....

.....

.....

.....

.....

.....

.....

.....

.....

.....

.....

.....

.....

.....

.....

.....

.....

(How does a potentiometer function?)

The students should first recognize from a model that a given voltage can be divided with a potentiometer (voltage divider) and that the dividing function is based on the law for the series connection of resistances $U_1/R_1 = U_2/R_2 = \dots = U_n/R_n$.

For a homogen wire is $I_1/R_1 = I_2/R_2 = \dots = I_n/R_n$.

The potentiometer must hereby not be under load.

Subsequently to this, the students should vividly experience the functioning of an actual potentiometer by working with it.

Notes on Set-Up and Procedure

In the first experiment, the lengths l can be freely chosen. The measured values for l and U can be best compared with each other, however, when the wire lengths are in the ratio of approximately 4:3:2:1.

In the second experiment it must be ensured that the preset power supply voltage does not exceed 5 V, because of the load-carrying ability of the lamp.

Observations and Measurement Results

Table 1

| $\frac{l}{m}$ | $\frac{U}{V}$ | $\frac{U}{l} \frac{V}{m}$ |
|---------------|---------------|---------------------------|
| 0.202 | 1.00 | 5.0 |
| 0.158 | 0.78 | 4.9 |
| 0.102 | 0.51 | 5.0 |
| 0.055 | 0.28 | 5.1 |

- (1) During the turning of the knob from the 0 mark up to the stop, the voltage increases from 0 V to a maximum value; during turning back, it decreases again back down to 0 V.
- (2) During the turning of the knob from the 0 mark up to the stop, the lamp starts to glow feebly, becomes continually brighter and attains its maximum brightness when the potentiometer knob reaches its stop; during turning back, the brightness of the lamp gets slowly less and less until the lamp goes out.

Evaluation

1. Refer to Fig. 3. The lengths of wire used l and the corresponding voltages U are directly proportional to each other, as the points lie on a straight line.
2. Refer to column 3 of Table 1.
 $U/l = \text{constant}$. It follows from this that:
 $U \sim l$ or $U_1/l_1 = U_2/l_2 = \dots = U_n/l_n$.
3. A potentiometer can be used to divide a given voltage into smaller voltages as required.
4. Potentiometers are used in electronic devices as components for the adjustment of loudness and brightness. The power supply also has a potentiometer for the adjustment of the output voltage.


Remarks

Potentiometers of low load have a carbon layer instead of a resistor wire.

You should also use the clarifying term voltage divider during your introduction to the potentiometer.

When the sliding contact of the potentiometer is connected to the end of a resistor path, then the potentiometer can be used as a component for changing resistance.

Fig. 3

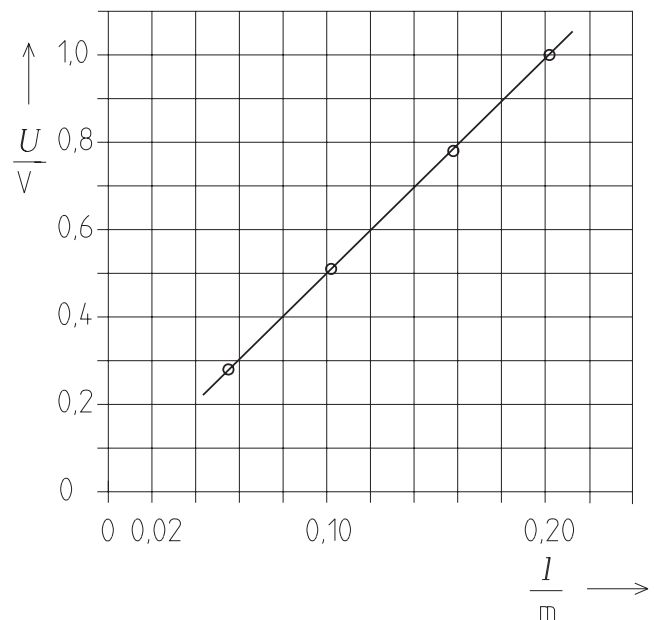

T**EEP
2.9**

The potentiometer



(How does a potentiometer function?)

Room for notes