

Steam pressure and atmospheric pressure (with support mate-rial)



Physics	Mechanics	Mechanics of liquids & gases	
Difficulty level	AA Group size	Preparation time	Execution time
easy	1	10 minutes	10 minutes

This content can also be found online at:



http://localhost:1337/c/5fd78c04ce3751000353f2d9





PHYWE



General information

Application PHYWE



Setup

The temperature dependent behaviour of fluids and gases has many implications. The increased pressure, that emerges when fluids turn to gases, can have dangerous consequences in enclosed spaces.



Other information (1/2)

PHYWE



Prior

knowledge



Main

principle

No prior knowlege is required.

Water is heated in a container until it boils, it is tightly sealed following this. The pressure in the container decreases as it cools. It is pressed together by the external air pressure.

Other information (2/2)

PHYWE



Learning

objective



Tasks

The goal of this experiment is to learn about the strong forces involved in the temperature dependent behaviour of fluids and gases.

1. Use the burner to test the Thermal expansion and thermal contraction in a tin can.



Equipment

Position	Material	Item No.	Quantity
1	Tripod base PHYWE	02002-55	1
2	Support rod, stainless steel, 750 mm	02033-00	1
3	Right angle clamp expert	02054-00	1
4	Support ring, i.d. 130mm,	37721-03	1
5	Tin can, 1 litres	02678-02	1
6	Beaker, 100 ml, plastic (PP)	36011-01	1
7	Wire gauze with ceramic, 160 x 160 mm	33287-01	1
8	Butane burner, Labogaz 206 type	32178-00	1
9	Butane cartridge C206, without valve, 190 g	47535-01	1





PHYWE



Setup and Procedure

Setup PHYWE

- Attach the support ring to the support rod using the right-angle clamp, place the wire gauze, set the height in such a way that the distance to the burner amounts to approx. 1 cm.
- Fill approx. 50 ml of water into the tin can and place this on the wire gauze.
- Place the screw top on the opening nozzle of the can but do not seal it yet.



Fig. 1





Procedure PHYWE

- Heat the tin can on a low flame.
- Allow the water to boil for a few minutes (the air should completely escape from the can if possible).
- Remove the burner and switch it off.
- Screw the screw top of the tin can tight (cloth, gloves).
- Let the tin can cool down.

Attention! Do not touch the hot tin can. Danger of getting burned!



Fig. 2

PHYWE



Evaluation





Results

When heating the water begins to boil then the steam escapes beneath the lid that has not been screwed tight. When it cools down the tin can is suddenly pressed together after a few minutes. This generally occurs in two or three steps.

Remarks:

To be able to estimate the pressure or forces that arise the vapor pressure of water at various temperatures must be known (table 1).

At an atmospheric pressure of 1013 hPa (normal pressure) water boils at 100°C, the vapor pressure at this temperature is thus the normal pressure. Once the steam cools to 20°C for instance the pressure in the canister is a mere 23 hPa. The atmospheric pressure is thus approx. 40 times greater than the vapor pressure in the vessel's interior.

The tin can has thin walls and additionally has edges on its lower and upper sides. It can therefore be dented in the event of this great pressure difference.

Evaluation PHYWE

If a liquid vaporizes or evaporates steam is present above its surface. This spreads out in the entire space that is available. The water boils for a few minutes so that there is no more air in the can (if possible) and instead just water and steam. The liquid vapor has a vapor pressure that is only dependent on the temperature. It is the same volume as the ambient air pressure when the liquid boils. The water and steam cools down again if the burner is removed. As a result the pressure (vapor pressure) in the interior of the canister decreases and the difference to the external air pressure is so great that the canister suddenly dents at one or several points.

T [°C]	[hPa]
20	23
100	1013

Table 1: Temperature T and vapor pressure p of water

