

PHYWE Systeme GmbH & Co. KG
Robert-Bosch-Breite 10
D-37079 Göttingen

Telefon +49 (0) 551 604-0
Fax +49 (0) 551 604-107
E-mail info@phywe.de
Internet www.phywe.com

Operating instructions

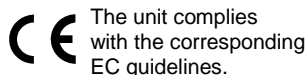


Fig. 1: Diodelaser, 1 mW, 532 nm (green) with short stem 08764-99

TABLE OF CONTENTS

- 1 SAFETY PRECAUTIONS
- 2 PURPOSE AND CHARACTERISTICS
- 3 DESCRIPTION
- 4 HANDLING
- 5 OPERATING INSTRUCTIONS
- 6 TECHNICAL DATA
- 7 WASTE DISPOSAL

1 SAFETY PRECAUTIONS



Warning of Laser beams

Laser radiation
Do not expose to the beam
Laser Class 2
(according to DIN EN 60825-1)

- Never look into the laser beam while the laser is switched on. Permanent eye damage may occur.

- Never direct the beam towards the face or eyes of persons, regardless of their distance.
- When using the laser, make sure that a functioning visual protection against radiation escaping from the test setup is guaranteed.
- Do not start up this instrument in case of visible signs of damage to it.
- Only use the instrument for the purpose for which it was designed.
- Only operate the laser with the supplied power supply unit.
- Never look directly into the laser beam, not even if it is expanded in its diameter, e.g. by upstream lenses. The optical phenomena to be examined should always be observed on a projection or transparent screen where the irradiance of the observing eye is reduced to a safe level by diffuse reflection or scattering. All optical devices in the beam path shall be mounted in such a way that an unintentional change of their position is not possible; no persons shall be endangered, e.g. by reflection of the beam, which can also occur on objects not belonging to the test setup.

2 PURPOSE AND CHARACTERISTICS

This diode-pumped, frequency-doubled YAG solid-state laser provides a compact monochromatic light source that is particularly suitable for experiments on interference and diffraction. Since the human eye is particularly sensitive to green light, the use of this laser is preferable to a red laser of equal power.

The laser meets the DIN safety requirements of laser protection class 2 and is therefore approved for schools.

3 DESCRIPTION

The lightproof housing used reliably prevents the escape of uncontrolled radiation. The connection to the mains is made via the plug-in power supply unit included in the scope of delivery.

To set up the laser in a tripod base or the rider of an optical bench, a 10 mm handle (included in the scope of delivery) is inserted into the housing.

4 HANDLING

Before starting the laser, hold it in a tripod foot or in the rider of an optical bench. The laser is to be set up in such a way that the laser beam cannot hit anyone in the eyes (see point 1, danger notes). The laser may only be connected to a mains socket using the supplied power supply (max. output voltage 3.3 VDC).

5 OPERATING INSTRUCTIONS

This high-quality instrument fulfils all of the technical requirements that are compiled in current EC guidelines. The characteristics of this product qualify it for the CE mark.

This instrument is only to be put into operation under specialist supervision in a controlled electromagnetic environment in research, educational and training facilities (schools, universities, institutes and laboratories).

This means that in such an environment, no mobile phones etc. are to be used in the immediate vicinity. The individual connecting leads are each not to be longer than 2 m.

The instrument can be so influenced by electrostatic charges and other electromagnetic phenomena that it no longer functions within the given technical specifications. The following measures reduce or do away with disturbances:

Avoid fitted carpets; ensure potential equalization; carry out experiments on a conductive, earthed surface, use screened cables, do not operate high-frequency emitters (radios, mobile phones) in the immediate vicinity.

6 TECHNICAL DATA

(typical for 25°C)

Operating temperature range: 15... 25

Laser


Laser type:	Diode
Wavelength:	532 nm not polarised
Power:	1 mW
Laser protection class:	Class 2
Beam diameter (at output):	0.5 mm
Beam divergence:	1.2 mrad
Connection for accessories (e.g. for beam expansion):	M20 external thread
Handle	
Length	75 mm
Diameter:	10 mm

Power supply unit

Input voltage range:	100...240 V~
Mains frequency	50/60 Hz
Output voltage	3.3 VDC
Output current	150 mA
DC-socket:	2.1 mm

7 WASTE DISPOSAL

The packaging consists predominately of environmentally compatible materials that can be passed on for disposal by the local recycling service.

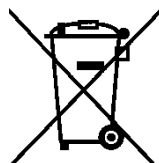

Should you no longer require this product, do not dispose of it with the household refuse.

Please return it to the address below for proper waste disposal.

PHYWE Systeme GmbH & Co. KG
Abteilung Kundendienst (Customer Service)
Robert-Bosch-Breite 10
D-37079 Göttingen

Phone +49 (0) 551 604-0
Fax +49 (0) 551 604-107