

PHYWE Systeme GmbH & Co. KG  
Robert-Bosch-Breite 10  
D-37079 Göttingen

Telefon +49 (0) 551 604-0  
Fax +49 (0) 551 604-107  
E-mail info@phywe.de  
Internet www.phywe.com

### Operating instructions

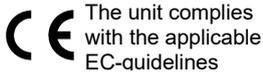


Fig. 1: 12951-00 Cobra SMARTsense Light & Color

## TABLE OF CONTENTS

- 1 SAFETY PRECAUTIONS
- 2 PURPOSE AND CHARACTERISTICS
- 3 FUNCTIONAL AND OPERATING ELEMENTS
- 4 NOTES ON OPERATION
- 5 HANDLING
- 6 TECHNICAL DATA
- 7 SCOPE OF DELIVERY
- 8 ACCESSORIES
- 9 CONFORMITY
- 10 DISPOSAL
- 11 NOTES ON BATTERY AND RECHARGEABLE BATTERY DISPOSAL

## 1 SAFETY PRECAUTIONS



### Caution!

- Carefully read these operating instructions completely before operating this instrument. This is necessary to avoid damage to it, as well as for user-safety.
- Only use the instrument for the purpose for which it was designed.
- Only use the instrument in dry rooms in which there is no risk of explosion.
- Protect the instrument from dust, moisture and vapours. Use a slightly moist lint-free cloth to clean the instrument. Do not use aggressive cleaning agents or solvents.
- Take care that no liquid penetrates in through the housing openings, as such penetration would result in damage to Sensor.
- Do not open the unit.
- Avoid pointing the white LED directly at your eye. The light is bright and may cause discomfort.

## 2 PURPOSE AND CHARACTERISTICS

The sensor is used to measure light in the visible and ultraviolet range, as well as to detect colour components and transmit the measured values via Bluetooth or USB to any terminal devices such as tablets, smartphones, etc.

### 3 FUNCTIONAL AND OPERATING ELEMENTS

#### 3.1 Operating elements

The sensor has an on-button and two LEDs for indicating the Bluetooth and battery charge status.

##### Power Button

Short pressed	Switch on white LED
Long pressed (>3s)	Switching the device on/off
Pressed 3x quickly	Start offline measurement
Pressed 2x quickly	Stop offline measurement

##### Bluetooth-LED

Flashing red every 2 seconds	Not connected
Flashing green every 2 seconds	Connected to the terminal device
Flashing green every 4 seconds	Running measurement

##### Battery charge LED

Flashing red every 2 seconds	Low battery
Illuminated red	Active charging process
Illuminated green	Charging process completed

#### 3.2 Functional elements

On the right side of the sensor are the sensor elements for light, UV and RGB/IR as well as a white illumination LED.

#### 3.3 USB port

The battery, which is permanently installed in the sensor, is charged via the type C USB port. Furthermore, communication with a computer takes place via this interface.

### 4 NOTES ON OPERATION

This device fulfils all of the technical requirements that are compiled in current EC guidelines. The characteristics of this product qualify it for the CE mark. The individual connecting leads are each not to be longer than 2 m.

The instrument can be so influenced by electrostatic charges and other electromagnetic phenomena (HF, bursts, indirect lightning discharges) that it no longer works within the given specifications. Carry out the following measures to reduce or eliminate the effect of such disturbance: Ensure potential equalization at the PC (especially with Laptops). Use screening. Do not operate high frequency emitters (e.g. radio equipment or mobile radiotelephones) in the immediate vicinity. When a total failure of the instrument occurs, unplug it and plug it back in again for a reset.

### 5 HANDLING

This section describes the start-up of the sensor and the recording of measurement data. Please read this section thoroughly in order to avoid failures or operating errors.

#### 5.1 Charging process

Use a USB-C cable to connect the sensor to a computer or USB charger (not included).

During the charging process, the battery charge LED lights up red. When the charging process is complete, the battery charge LED lights up green.

The charging time for a completely discharged battery is 3 hours maximum.



Disconnect the charger at the latest four hours after the completion of the charging process. Otherwise, the service life of the battery may be negatively affected.

#### 5.2 Start-up

Switch on the sensor by pressing the power button for more than 3s. Now the Bluetooth LED flashes red. Start the software and select the sensor.

If the sensor is to be used via the USB interface, it does not need to be switched on. The sensor is connected directly to the end device using the supplied USB cable.

There is a 9-digit code on the back of the sensor (Fig.2). The last 4 digits of the code are displayed as the sensor name in the software (Fig.3). This enables the precise assignment of the sensors within the software.



Fig. 2



Fig. 3

#### Selection of the sensor via the Bluetooth interface

Make sure that the Bluetooth interface is activated on the terminal device (PC/Tablet/Smartphone) and that the software is allowed to access the interface.

After the sensor has been selected in the software, the LED flashes green to indicate that the connection has been established correctly. After the sensor has been coupled with the software, the sensor is no longer visible to other users in the software, and therefore can no longer be selected.

If the sensor is switched on and not connected, it switches off automatically after 5 minutes.

#### Selection of the sensor via the USB interface

For this purpose the sensor must be plugged into the USB port of the end device. It is not necessary to switch on the sensor. The sensor is automatically recognized and displayed. It can be selected and connected directly.

#### 5.3 Offline measurement

Switch on the sensor by pressing the power button for more than 3s. To start an offline measurement, press the power button 3 times in quick succession. The Bluetooth LED then flashes green 3 times in rapid succession to acknowledge the successful start. To stop a measurement, press the switch-on button 2x in quick succession. The Bluetooth LED also acknowledges this by flashing quickly.

Offline measurements can be read out via the measureAPP or measureLAB software. Furthermore, offline parameters such as data rate and measurement duration can be set. After the set measurement duration has elapsed, the offline measurement is automatically terminated. However, the measurement can always be ended prematurely by pressing the switch-on button.

## 5.4 Switching the white illumination LED on/off

Press the switch-on button briefly to switch on the LED. Pressing the switch-on button again switches the LED off.



Avoid pointing the white LED directly at your eye. The light is bright and may cause discomfort.

## 6 TECHNICAL DATA

Operating temperature range: 5 - 40°C  
Rel. humidity < 80%

### Illuminance

Measuring range	0 ... 200 kLx
Resolution	1 Lx
Max. data transfer rate	1000 Hz

### UV radiation (UVA, UVB, UVC)

Measuring range	0 ... 400 W/m <sup>2</sup>
Resolution	0,1 W/m <sup>2</sup>
Wavelength	210 ... 370 nm
Max. data transfer rate	1 Hz

### RGB/IR

Measuring range	0 ... 1000 (rel)
Wavelength	
- Red	580-720 nm
- Green	450-600 nm
- Blue	400-520 nm
- Infrared (IR)	780-1100 nm
Max. data transfer rate	1 Hz

Max. wireless range (open field)	30 m
Battery	LiPo 3.7 V / 2000 mAh
Dimensions (w x h x d)	92 x 37 x 25 mm
Weight	60 g

## 7 SCOPE OF DELIVERY

The extent of delivery is as follows

- Cobra SMARTsense Light & Color 12951-00
- USB connecting cable type C 07935-00
- Tripod rod with thread (10cm)
- Operating instructions

## 8 ACCESSORIES

- USB connecting cable type C 07935-00
- Cobra SMARTlink 12998-99
- USB-Bluetooth-Adapter 07936-00
- Software measureLAB 14580-61
- Free measureApp available from supplier portals

iOS

Android

Windows



## 9 CONFORMITY



PHYWE Systeme GmbH & Co.KG hereby declares that the radio system type 12951-00 complies with the 2014/53/EU directive. The complete text of the EC Declaration of Conformity is available at the following Internet address:

[www.phywe.com/en/ec-declaration](http://www.phywe.com/en/ec-declaration)

## 10 DISPOSAL

The packaging mainly consists of environmentally-friendly materials that should be returned to the local recycling stations.



Do not dispose of this product with normal household waste. If this unit needs to be disposed of, please return it to the address that is stated below for proper disposal

PHYWE Systeme GmbH & Co. KG  
Customer service department  
Robert-Bosch-Breite 10  
D-37079 Göttingen

Phone +49 (0) 551 604-0  
Fax +49 (0) 551 604-107

## 11 NOTES ON BATTERY AND RECHARGEABLE BATTERY DISPOSAL

As we sell batteries and rechargeable batteries or devices containing batteries and rechargeable batteries, we are obliged under the Battery Act to inform you of the following: Batteries and rechargeable batteries may not be disposed of with household waste, but you are legally obliged to return used batteries and rechargeable batteries. Used batteries may contain harmful substances that can damage the environment or your health if they are not stored or disposed of properly. Batteries also contain important raw materials such as iron, zinc, manganese or nickel and are recycled. You can either send the batteries back to us after use or return them free of charge in the immediate vicinity (e.g. in shops or at municipal collection centres). Batteries or rechargeable batteries that contain harmful substances are labelled with the symbol of a crossed-out dustbin.