HORO WÄRMESCHRÄNKE

HORO Dr. Hofmann GmbH, Rudolf- Diesel- Str. 8, 73760 Ostfildern
Tel ++ 49 / 711 / 341 10 59 Fax ++ 49 / 711 / 348 27 08 info@horo-dr-hofmann.de

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Instructions for the HORO Electric Burner

Check that the connectors are of the correct voltage. Switch on with the toggle switch at the connector. After about 2 minutes the furnace will be glow red. Full working capacity can only be reached when connected to the correct voltage (voltage drops will decrease the temperature).

Burner components such as the white porcelain wall or the soldering iron holder can reach temperatures in excess of 60°C (even after switching off).

Burners that are damaged are not to be used. The burner is to be set up as far away as possible from combustible materials.

Avoid covering of the burner to facilitate adequate ventilation to the furnace area.

Children do not recognize the dangers of electrical heating devices and should therefore be permitted use only under direct adult supervision.

Repairs are to be undertaken by the manufacturer and only with original replacement parts. We will assume no guarantee or responsibility for unauthorized opening or repair of the burner.

All burners are manufactured to EEC guidelines.

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Hereby we declare, that this product is in conformity with the following standards or normative documents with respect to the design:

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¹ Der GmbH- Geschäftsführer ist für Zusammenstellung der technischen Unterlagen bevollmächtigt.
The CEO of the GmbH is authorized for publishing the technical documentation